

## REFERENCE ELECTRODE INSTRUCTIONS

---

**DO NOT TWIST CAP!**

**Push  
Middle  
Section  
DOWN to  
Open the  
Fill Hole**



**DO NOT TWIST CAP!**

---

Upon receipt from the factory, the reference electrode contains a small amount of distilled water. It is not necessary to remove this water before filling the electrode.

Before using the electrode, fill the inner chamber with the Fill Solution (provided). After filling, place the electrode in a small beaker of water (preferably distilled). This will permit the Fill Solution to flow out through the bottom of the electrode and establish an ionic conduction path through the sintered glass frit at the bottom of the electrode.

Reference Fill Solution  
is available from  
**Analytical Sensors, Inc.**  
1-800-545-6132  
[www.asi-sensors.com](http://www.asi-sensors.com)